

Latest development of α emitter imaging and quantification on a large Field Of View

Arthur Bongrand, Samuel Duval, Jérôme Donnard, Julie Champion

▶ To cite this version:

Arthur Bongrand, Samuel Duval, Jérôme Donnard, Julie Champion. Latest development of α emitter imaging and quantification on a large Field Of View. International Symposium on Trends in Radiopharmaceuticals (ISTR-2023), Apr 2023, Wien, Austria. hal-04098723

HAL Id: hal-04098723 https://hal.science/hal-04098723

Submitted on 16 May 2023

HAL is a multi-disciplinary open access archive for the deposit and dissemination of scientific research documents, whether they are published or not. The documents may come from teaching and research institutions in France or abroad, or from public or private research centers. L'archive ouverte pluridisciplinaire **HAL**, est destinée au dépôt et à la diffusion de documents scientifiques de niveau recherche, publiés ou non, émanant des établissements d'enseignement et de recherche français ou étrangers, des laboratoires publics ou privés.

Latest development of a emitter imaging and quantification on a large Field Of View



1: Laboratoire Subatech, IMT Atlantique, Nantes Université, CNRS/IN2P3, France

2: S.A.S. Ai4R, France

arthur.bongrand@subatech.in2p3.fr



Université



#273

Currently, the characterization of the nature and spatial distribution of radionuclides in a sample is time consuming and fastidious. Indeed, it necessarily requires the use of two distinct analytical techniques and detectors of different nature.

Moreover, the spatial resolutions of the two methods are often detector dependent, which makes it difficult to integrate and interpret these two measurements together.

Can we perform a fast and accurate quality control of radiopharmaceuticals with a single system?



Digital Autoradiography...

AEA ISTR

International Atomic Energy Agency

\rightarrow an accurate up-to-date emission imaging technique !

- Allows to study and locate in 2 dimensions the radioactivity emanating from a surface sample at scales until **10 micrometer.**
- Well suited for pre-clinical TAT as it allows understanding and quantifying biodistributions \checkmark in tissues at high spatial resolution and low activity.
- Digital autoradiography \rightarrow Ability to analyze and reanalyze endlessly the digital data acquired!
- Variety of applications: Radio-TLC, Geological Studies, Radio Toxicology...

#275

Time, t/T1/2 of daughter

BeaQuant: a particle by particle digital autoradiography system





 \rightarrow Usually the quality controls are made from a 1D projection where we can distinguish the free ²²⁵Ac which remains on the deposit spot from the labelled ²²⁵Ac which migrates. To limit interpretation errors, the analysis must be performed at **secular equilibrium** (~7h).



→ The STSAA reduces the time needed for a quality control after elution to only one hour allowing both isotope localization and identification. Thus, the purity is deduced from the ratio of

The Spatio Temporal Spectroscopic Autoradiography Analysis (STSAA)



a radioactive nuclide decays at the same rate as it is produced.

- \rightarrow The concentration of daughter nuclei in the radioactive equilibrium can be modeled at any time using the **Bateman equation** provided that the characteristics of the decay chain and if the initial concentrations are known.
- \rightarrow For our application, as these initial concentrations are unknown, we have developed a method to calculate these concentrations from the evolution of the measured count rate.



free to labeled ²²⁵Ac.

Discussion

Digital autoradiography combined with a **temporal spectroscopy** method allows the **quick** and accurate identification of different radioisotopes and to observe their localization.

The increased robustness of this method comes from its low dependence on the operator's choices compared to 1D method.

From the results, we can observe the **contribution** of a radioisotope called "free ²¹³Bi". Clearly visible on the migration front, free ²¹³Bi corresponds to ²¹³Bi which is not radiolabeled and whose parent, ²¹⁷At, has not decayed locally. The most probable hypothesis is that the free ²¹³Bi comes from the mother solution. This free ²¹³Bi has also been observed in biological samples. Studies are underway to understand its biological impact.

Mouse kidney labelled with ²²⁵Ac-PSMA

≥ 0.14

⊈ 0.12

0.1

<u>9</u> 0.08

는 0.06

ပ<u>ိ</u> 0.04

0.02





Conclusions

Digital Autoradiography with spectroscopic method:

Promising in the \checkmark for nuclear medicine applications like quality control of radionuclides of radiopharmaceuticals !

- A precise tool to understand the temporal behavior of a chain of labeled \checkmark isotopes.
- Precise image & quantification from mBq to kBq.

References

- -

[1] H. Lefeuvre et al., « Spectroscopic autoradiography of alpha particles using a parallel ionization multiplier gaseous detector », Nuclear Instruments and Methods in Physics Research Section A: Accelerators, Spectrometers, Detectors and Associated Equipment, vol. 1035, p. 166807, 2022.

[2] R. Eychenne et al., « Overview of the Most Promising Radionuclides for Targeted Alpha Therapy: The "Hopeful Eight" », Pharmaceutics, vol. 13, n°6, Art. no 6, 2021.

Acknowledgments

We thank the GIP Arronax for providing of their facilities and a ²²⁵Ac-PSMA source, especially Dr. Cyrille Alliot for his excellent expertise in the field of radiopharmaceutical quality control.

More experiments !

Perspectives

acquisition of nuclear data for the production \rightarrow Perform simultaneous comparisons with standard quality control methods.

-> Perform multiple quality controls of the same batch simultaneously using the BeaQuant's large Field Of View. Tb 149

 \rightarrow Explore the method for other radiopharmaceuticals [2]

At 211 4.1h 7.2h α 3.<mark>97</mark> β⁺ γ 352; 165.



International Symposium on Trends in Radiopharmaceuticals (ISTR-2023), CN-310 Vienna, Austria; 17-21 April 2023

